

AN033

Adaptive Optics Course Outline

Author: Justin Mansell, Ph.D.

Revision: 9/9/11

Introduction

We have sold adaptive optics (AO) kits to many universities. Our AO hardware and software is perfect for students learning about AO because it is easy to get something simple working immediately and is also flexible enough for students to write any part of the AO control system themselves using our provided Matlab class interface. In this application note we provide an example outline for teaching a graduate course in adaptive optics either using AOS hardware or a modeling-only based approach.

Outline for a Course in

Adaptive Optics with a Laboratory

Requirements: Electromagnetics Physics, Calculus, Introductory Computer Programming

Recommended Prerequisite Course(s): Fourier Optics, Introductory Optics

Recommended Reference Texts: J. Goodman's Fourier Optics, A. E. Siegman's Lasers, J. Schmitt's Optical

Modeling, Application Notes at www.aos-llc.com

Recommended Textbook: R. Tyson's Introduction to Adaptive Optics

Week	Lecture	Laboratory
1	Introduction to Adaptive Optics	Introduction to Matlab
	Applications of Adaptive Optics	
2	Review of Maxwell's Equations	Introduction to Matlab Script & Function
	Derivation of the Wave Equation	Programming
3	Review/Introduction to Fourier Optics	Write a script to perform 1D Fourier
		Transforms in Matlab
4	Fresnel & Fraunhofer Diffraction	Setup a large collimated beam using a shear
	Gaussian Beam Propagation	plate.
	ABCD Propagation	Measure Gaussian Beam Propagation in the
		laboratory with a beam profiler or a camera.
		Alternative Laboratory:
		Write a Matlab script to model the
		propagation of a Gaussian beam through a

		series of lenses using ABCD matrix theory.
5	Numerical Methods for Modeling Optical Propagation Fourier Mesh Requirements	Write a function to propagate an arbitrary 2D optical field.
6	Siedel Polynomials & Zernike Polynomials Effect of Aberrations on a top-hat beam propagating to focus.	Tilt a lens and measure the induced aberration on the focus of a beam. Place a beam splitter in the converging beam. Tilt the beam and show the induced aberration. Model the propagation of a beam with an aberration in Matlab. Compare the lab results with your model
		Alternative Lab: Write a function for the real Zernike polynomials. Show the effect of each of the first 15 polynomials on the far-field of a circular top-hat beam.
7	Wavefront Sensors – Interferometers, Hartmann Sensors, and Curvature Sensors	Setup a collimated beam into a Hartmann wavefront sensor. Calibrate the wavefront sensor on the beam. Measure a lens placed into the beam. Observe the effect of moving the lens away from the wavefront sensor. Observe the effect of tilting the lens. Setup a simple 2-lens telescope and observe the effect of moving the lenses. Alternative/Additional Lab: Build an interferometer using a HeNe laser, a beam splitter or sampler, and 2 mirrors (one of which could be deformable). Alternative Lab: Model the effect of a Zernike aberration on a variety of wavefront sensors (interferometer,
8	Deformable Mirrors – Plate-type, Membrane, Bimorph, & MEMS Comparison of the influence functions Comparison of the advantages and disadvantages for common AO applications.	Hartmann sensor, curvature sensor) Reflect a collimated beam from the membrane deformable mirror. Observe the effect of electrically adjusting the actuators on the beam propagated a distance from the DM. Observe the effect of the actuators near the focus of a beam. Alternative Lab: Perform the experiment above numerically

		using a simulated DM.
		Alternative Lab 2:
		Choose an application of adaptive optics
		(ophthalmology, directed energy, imaging,
		astronomy) and make an argument for which
		DM would be best for this application now
		and in the future.
9	Kolmogorov Turbulence	Measure turbulence in the lab by
	Statistics of Kolmogorov Turbulence	propagating a collimated beam over a large a
	Phase Structure Functions	distance as is practical and observe the effect
	Create a Matlab function to generate	of the air motion.
	Kolmogorov turbulence using the	Place an arm under the beam and observe
	prophability density function.	the air motion due to the heat coming off of
	Greenwood frequency	the arm.
	Statistics of compensation of Kolmogorov	Place a warm/hot object (coffee cup with
	turbulence using a DM	warm water, hair dryer, hotplate, etc.)
	tarbalence asing a bivi	underneath the beam and measure the
		induced aberration again.
		Use Matlab to analyze the data and calculate
		the phase structure function for each case
		the phase structure function for each case
		Alternative Lab:
		Use Matlab to model generation and
		compensation of Kolmogorov turbulence
		with a DM with Gaussian influence functions.
10	Metric Adaptive Optics	Use camera or photodiode feedback to close
10	Common Metrics: Sharpness, Second	a metric AO loop to minimize spot size
	Moment, RMS Slopes, Beam Shape, etc.	(second moment).
	Common Techniques: Stochastic Parallel	(Second Moment).
	Gradient Descent (SPGD), Guided	Alternative Laboratory:
	Evolutionary Simulated Annealing (GESA),	Take a provided aberrated optical field and
	Axial Searching, Stochastic Axial Searching	use metric AO techniques to compensate the
	Total Scaroming, Scooliastic Total Scaroming	beam with either a set of Zernike
		polynomials or with DM influence functions.
11	Matrix-Based AO	Add a wavefront sensor to the experiment
	Simple tilt control	from the prior week. Perform active
	SVD Inverse with Mode Removal	flattening of a membrane DM. Compare the
	Slope Discrepancy	results with the metric AO result.
	Ciopo Dissi oparis,	7.55.8.65 17.67 18.77 18.
		Alternative Laboratory:
		Take the same field from the prior week and
		compensate it using a modeled wavefront
		sensor feedback.
12	Advanced Concepts / State of the Art	Advanced Concepts / State of the Art
13	Review for the Final Exam or	Final Exam or Research Paper or Final Oral
15	More Advanced Concepts	Presentation (15 minutes each)
	iviore Advanced Concepts	r resentation (15 minutes each)